

# Technical Data Sheet

Doc No Ab/TDS/Eqpt-241

Rev No : 05 Date 02/2022-23

## TECHNICAL DATA SHEET

Abdos Labtech Pvt. Ltd.

### WAVEX – CLASSIC TUBE ROTATOR

Valid for Catalog No. E11300

#### 1. Description/Specification

##### 1.1 Description:

1.1.1 Wavex tube rotators provide gentle but effective mixing of biological samples, held in 1.5 ml to 50 ml tubes.

1.1.2 Used in a variety of applications including immune precipitations, prevention of blood coagulation & latex diagnostics.

1.1.3 Wavex-Classic Tube Rotator has Disk Type Design.

##### 1.2 Design:

Disk type design

##### 1.3 performance Specifications:

1.3.1 Voltage (AC)	: 220~240 Volt
1.3.2 Frequency (Hz)	: 50
1.3.3 Power (W)	: 20
1.3.4 Tilt Angle	: 0- 90°
1.3.5 Motor type	: DC Motor
1.3.6 Speed range (RPM)	: 0 ~ 80
1.3.7 Operation time	: Continuous
1.3.8 Dimensions	: 300x220x310mm
1.3.9 Permissible ambient Temp	: 5- 40°
1.4.0 Permissible Relative humidity:	80 %
1.4.1 Speed Display :	Scale

##### 1.4 Material: Body – ABS, Stainless Steel



Abdos Labtech Private Limited  
(A Division of Abdos Group)  
(An ISO 9001,13485,14001,18385,GMP & CE certified company)  
Shiv Ganga Industrial Estate  
Plot No-1, Lakheswari, Bhagwanpur,  
Roorkee-247667, (UK) INDIA

Registered office  
13A Government  
Place East, 2nd Floor Kolkata  
West Bengal – 700069  
CIN: U85100WB2008PTC123903  
Corporate Office  
319, 3rd Floor, DLF Tower A, DDA Dist. Centre  
Jasola, New Delhi - 110025

Ph. No: +91- (0) 1332235277  
Email:labtech@abdosindia.com  
Web: [www.abdoslivesciences.com](http://www.abdoslivesciences.com)  
GST: 05AAGCA8605A1Z7

# Technical Data Sheet

**1.5 Color:** Dark Blue

**1.6. Spare Accessories:**

E11301 Plastic clips for 1.5ml x 12 MCT, Disk accessory – 12pcs/case

E11302 Plastic clips for 15ml x 12 MCT, Disk accessory – 12pcs/case

E11303 Plastic clips for 50ml x 12 MCT, Disk accessory – 12pcs/case

1.7. CE certified

1.8 **Chemical resistance:** see website ([www.abdoslifesciences.com](http://www.abdoslifesciences.com))

**Warranty:** 12 months from the date of Installation or 15 months from the date of Invoicing

## Instruction for Use

Storage temp. Max 30°C

Before using the product check crackness, etc.

Check all the information printed on pouches or boxes.

## WARNING

**DO NOT** store strong oxidizing agents.

**DO NOT** place any plastic labware in a flame.

**DO NOT** mix any chemicals that may result in a thermic reaction, which can cause product failure

Consult local fire codes prior to storage of flammable liquids.



**Abdos Labtech Pvt. Ltd.**

Plot No. 1 Lakeshwari Bhagwanpur Roorkee 247667 (UK) India  
Email : [labtech@abdosindia.com](mailto:labtech@abdosindia.com), [www.abdoslifesciences.com](http://www.abdoslifesciences.com)  
Ph: +(91) 11 41081897



Electronic IFU



Abdos Labtech Private Limited  
(A Division of Abdos Group)  
(An ISO 9001,13485,14001,18385,GMP & CE certified company)  
Shiv Ganga Industrial Estate  
Plot No-1, Lakhewari, Bhagwanpur,  
Roorkee-247667, (UK) INDIA

**Registered office**  
13A Government  
Place East, 2nd Floor Kolkata  
West Bengal – 700069  
CIN: U85100WB2008PTC123903  
**Corporate Office**  
319, 3rd Floor, DLF Tower A, DDA Dist. Centre  
Jasola, New Delhi - 110025

Ph. No: +91- (0) 1332235277  
Email:[labtech@abdosindia.com](mailto:labtech@abdosindia.com)  
Web: [www.abdoslifesciences.com](http://www.abdoslifesciences.com)  
GST: 05AAGCA8605A1Z7

# Technical Data Sheet



**Abdos Labtech Private Limited**  
(A Division of Abdos Group)  
(An ISO 9001,13485,14001,18385,GMP & CE certified company)  
Shiv Ganga Industrial Estate  
Plot No-1, Lakheswari, Bhagwanpur,  
Roorkee-247667, (UK) INDIA

**Registered office**  
13A Government  
Place East, 2nd Floor Kolkata  
West Bengal – 700069  
CIN: U85100WB2008PTC123903

**Corporate Office**  
319, 3rd Floor, DLF Tower A, DDA Dist. Centre  
Jasola, New Delhi - 110025

**Ph. No: +91- (0) 1332235277**  
**Email:labtech@abdosindia.com**  
**Web: [www.abdoslivesciences.com](http://www.abdoslivesciences.com)**  
**GST: 05AAGCA8605A1Z7**